

AH401B

4-Channel Picoammeter



User's Manual



All Rights Reserved
© CAENels d.o.o.
Rev. 1.2 – November 2010

This product is licensed by Sincrotrone Trieste S.C.p.A.



CAENels d.o.o.
Kraška ulica, 2
6210 Sežana – Slovenija
Mail: info@caenels.com
Web: www.caenels.com

Table Of Contents

1. INTRODUCTION.....	7
1.1 THE AH401B PICOAMMETER	7
1.2 THE AH401B AT A GLANCE	8
1.3 FEATURES	8
1.4 DATA FORMAT.....	10
1.5 OFFSET CALIBRATION.....	11
2. SOFTWARE COMMANDS	13
2.1 COMMAND SYNTAX	13
2.2 COMMAND LIST	15
2.2.1 “ACQ” Command.....	15
2.2.2 “BDR” Command.....	17
2.2.3 “BIN” Command.....	18
2.2.4 “GET” and “?” Command.....	19
2.2.5 “HLF” Command.....	20
2.2.6 “ITM” Command.....	22
2.2.7 “NAQ” Command.....	23
2.2.8 “RNG” Command.....	24
2.2.9 “SUM” Command	25
2.2.10 “TRG” Command.....	26
2.2.11 “VER” Command	28
2.3 COMMAND TABLE SUMMARY	29
3. COMMUNICATION MODULES	31
3.1 MODULE HARDWARE INSTALLATION	31
3.2 “ETHERPIGGY” MODULE.....	32
3.2.1 “ETHERPiggy” IP Address Assign.....	32
3.2.2 “ETHERPiggy” Software Configuration	35
4. I/O CONNECTORS.....	39
4.1 POWER CONNECTOR	39
4.2 I/O INTERFACE CONNECTOR	39
4.2.1 The CONV Signal.....	40
4.2.2 The TRIGGER/GATE Signal.....	40
5. TECHNICAL SPECIFICATIONS	41

Document Revision	Date	Comment
1.2	November 19 th 2010	Added NAQ Command Added SUM Command Modified TRG Command
1.1	August 6 th 2009	Corrected Binary Output Format
1.0	October 27 th 2008	First Release



Safety information - Warnings

Read over the instruction manual carefully before using the instrument.
The following precautions should be strictly observed before using the AH401B:

WARNING

- Do not use this product in any manner not specified by the manufacturer. The protective features of this product may be impaired if it is used in a manner not specified in this manual.
- Do not use the device if it is damaged. Before you use the device, inspect the instrument for possible cracks or breaks before each use.
- Do not operate the device around explosives gas, vapor or dust.
- Always use the device with the cables provided.
- Turn off the device before establishing any connection.
- Do not operate the device with the cover removed or loosened.
- Do not install substitute parts or perform any unauthorized modification to the product.
- Return the product to the manufacturer for service and repair to ensure that safety features are maintained

CAUTION

- This instrument is designed for indoor use and in area with low condensation.



The following table shows the general environmental requirements for correct operation of the instrument:

Environmental Conditions	Requirements
Operating Temperature	5°C to 45°C
Operating Humidity	20% to 85% RH (non-condensing)
Storage Temperature	-20°C to 70°C
Storage Humidity	5% to 90% RH (non-condensing)

1. Introduction

This chapter describes the general characteristics and main features of the AH401B picoammeter.

1.1 The AH401B Picoammeter

The AH401B is a 4-channel, 20-bit resolution, low noise, high performance picoammeter. It is composed of a particular charge-integration input stage for low current sensing, coupled with a 20-bit sigma-delta ADC converter including a noise reduction digital filter.

The AH401B performs current measurements from 50pA (with a resolution of 50 aA) up to 1.8 μ A (with a resolution of 1.8 pA), with integration times ranging from 1ms up to 1s. Moreover, each input channel has two parallel integrator stages, so that the current-to-voltage conversion can be performed continuously also during the ADC conversion, avoiding any dead time in the data throughput.

It also performs digitization of the acquired current data, thus strongly minimizing the transmission length of analog current signals from the detector and providing directly digital data output, which is easy to transmit over long distances without any noise pick-up.

The AH401B is housed in a light, practical and extremely compact metallic box that can be placed as close as possible to the current source (detector) in order to reduce cable lengths and minimize possible noise pick-up during the propagation of very low intensity analog signals. It is specially suited for applications where multi-channel simultaneous acquisitions are required, a typical application being the currents readout from 4-quadrant photodiodes used to monitor X-ray beam displacements. Low temperature drifts, good linearity and the very low intrinsic noise of the AH401B allow obtaining very high precision current measurements.

The AH401B with its modular communication capability allows the user to freely select the type of communication interface, allowing control of the instrument with different types of programming languages and/or operating systems. Currently available communication modules are: RS232, RS422/485, USB and Ethernet (TCP/IP and UDP).

The efficient integration realized in the AH401B of the whole acquisition, counting and digitization chain in a single compact element allows simplifying and streamlining to a great extent the task to read out at high speed and with extremely low noise X-ray detectors, while also greatly enhancing the price/performance ratio and its user-friendliness.



1.2 The AH401B at a Glance

In Fig. 1 the AH401B and its I/O connections can be easily seen:



Fig. 1 - The AH401B I/O Connections

1.3 Features

The AH401B input stage is based on a typical charge integrator topology, thus the instrument really measures the charge integrated during the selected integration time. Eight full scale capacitor ranges are available as reported in the following table:

	Full Scale
RNG 0	1.8nC
RNG 1	50pC
RNG 2	100pC
RNG 3	150pC
RNG 4	200pC
RNG 5	250pC
RNG 6	300pC
RNG 7	350pC

Whereas the integration time can be varied between 1ms and 1s in incremental steps of 100 μ s.

In order to convert the input data from “integrated charge” to current the following formula must be applied:

$$I_{IN} = \frac{Q_{IN}}{t_{INT}}$$

In the following table the maximum and minimum currents that can be measured with the AH401B are reported, along with the LSB resolution, for each capacitor range:

	Full Scale	Equivalent Current FS @1ms t_{INT}	LSB Resolution @1ms t_{INT}	Equivalent Current FS @1s t_{INT}	LSB Resolution @1s t_{INT}
RNG 0	1.8nC	1.8 μ A	1.8pA	1.8nA	1.8fA
RNG 1	50pC	50nA	50fA	50pA	50aA
RNG 2	100pC	100nA	100fA	100pA	100aA
RNG 3	150pC	150nA	150fA	150pA	150aA
RNG 4	200pC	200nA	200fA	200pA	200aA
RNG 5	250pC	250nA	250fA	250pA	250aA
RNG 6	300pC	300nA	300fA	300pA	300aA
RNG 7	350pC	350nA	350fA	350pA	350aA

A host PC is necessary in order to operate the AH401B and properly set the desired parameters (i.e. range and integration time) and to acquire the converted data. Please refer to the “Software Commands” chapter for a complete description of available commands, their purposes and syntax.

A fully developed LabView™ GUI is available, allowing to plot the AH401B outputs up to 500 Hz and performing all necessary operations to convert the current data into positional information for all most commonly used Beam Position Monitors geometry. It also performs calculation of statistical data on the sampled buffers of data (average and rms) and, most importantly, provides the user with a FFT function that is of uttermost importance when using the AH401B as a spectrum analyzer to characterize the vibrational properties of the X-ray beam. The FFT can extend up to 250 Hz, fully covering the critical 0-100 Hz range where most of the beamlines require extreme stability in order not to spoil the science data acquisition.

1.4 Data Format

Since the AH401B works in “charge integrator mode” the data output values depend on the range and integration time selected. For this reason it is important to properly set these parameters in order to avoid input stage saturation and the corresponding data clipping (to 0 or to 1048575, the latter being the upper limit before saturation for the AH401B in raw digital data) in the output values.

Moreover, as an output offset has been built into the instrument to allow for the measurements of input signals near and below zero, this data offset must be taken into account for proper data conversion. For this reason an offset calibration procedure must be performed before using the AH401B. In the following table the correspondence between the read value and the full scale range is reported:

Data Value	Input Signal
1048575	Full Scale
4096	0
0	-0.4% Full Scale

The data read from the AH401B are the raw values of the 20-bit ADC converter, thus the data conversion from the read values to the corresponding current values must be performed by the user applying the following formula:

$$I_{IN} = \frac{FSR}{2^{20} - 1} \cdot (Value - Offset) \cdot t_{INT}$$

with:

- I_{IN} : the calculated input current;
- FSR: the selected Full Scale Range;
- Value: the raw data value from the AH401B;
- Offset: the data offset (~4096). For a proper estimation of this offset a calibration procedure must be performed (see “AH401B Offset Calibration” paragraph);
- t_{INT} : the selected Integration Time.

1.5 Offset Calibration

As explained in the previous paragraph, an output offset is built into the instrument to allow for the measurements of input signals near and below zero. For this reason with an input signal of 0 the read value from the AH401B is equal to ~4096. If a precise estimation of this offset is necessary, the following procedure must be followed:

- Place on each input connector a shielding cap in order to avoid noise pick-up (be aware that the central input pin must be left open and shielded but not grounded otherwise the read values are meaningless);
- Switch on the AH401B;
- Let it warm-up for about half an hour;
- Acquire a set of data (for example 100pts.) from each input channel and calculate the mean value;

The calculated mean values are the new Data Offset that must be taken into account in the current evaluation formula (see 1.4) that has to be applied to each channel.

1




2. Software Commands

This chapter describes the software commands used to tune and read the AH401B picoammeter.

2.1 Command Syntax

The command syntax used by the AH401B protocol is described in the following paragraphs.

The commands are sent in ASCII format and are composed by two fields, the “command field” and the “parameter field”, separated by a space. Instructions are not case sensitive and therefore the command string can freely use either uppercase or lowercase. Each instruction is terminated with a “carriage return” character (`\r` or 0x0D in hexadecimal notation), e.g.:

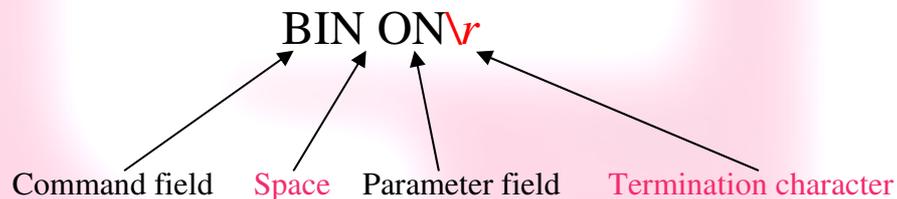

There are two types of software commands: “configuration commands” and “query commands”. The “query commands” usually have the same “command field” as the “configuration commands” with a “?” in the “parameter field”.

After each received command the AH401B replies with a string whose content depends on the type of command sent and its correctness.

There are three possible scenarios (with some exceptions, as described in the corresponding command paragraph):

1) *the command syntax is incorrect*: the AH401B will always answer with the string “NAK” terminated with a “carriage return/line feed” characters (`\r\n` or 0x0A0x0D in hexadecimal notation), e.g.:

`BIX ON\r` → ← `NAK\r\n`

example 2:

`BIN OOG\r` → ← `NAK\r\n`

2) *the command syntax is correct and it is a “configuration command”*: the AH401B will answer with the string “ACK” terminated with a “carriage return/line feed” (`\r\n`) characters, e.g.:

`BIN ON\r` → ← `ACK\r\n`

example 2:

`ITM 1000\r` → ← `ACK\r\n`

3) *the command syntax is correct and it is a “query command”*: the AH401B will answer with the echo of the “command field” and the value of the requested parameter, separated with a space and terminated with a “carriage return/line feed” (`\r\n`) characters, e.g.:

`BIN ?\r` → ← `BIN ON\r\n`

example 2:

`ITM ?\r` → ← `ITM 1000\r\n`


2.2 Command List

Please find hereafter the command list used by the AH401B with their syntax, the description of their purpose and the description of any special requirements.

2.2.1 “ACQ” Command


The purpose of the ACQ command is to start/stop the acquisition from the AH401B. As soon as the command is received the instrument starts to acquire data and immediately sends the generated digital data stream to the host PC. At power-up the instrument boots with the acquisition stopped (“ACQ OFF” - *default*), the user is thus required to start the data acquisition sending the “ACQ ON” command.



The type of data transmitted depends on the setting of the “binary mode” that can be changed using the “BIN” command (see “BIN Command” paragraph). With “BIN OFF” (*default*), data is sent in ASCII mode with the following format: “ch1 ch2 ch3 ch4\r\n”, with a space between each channel value:

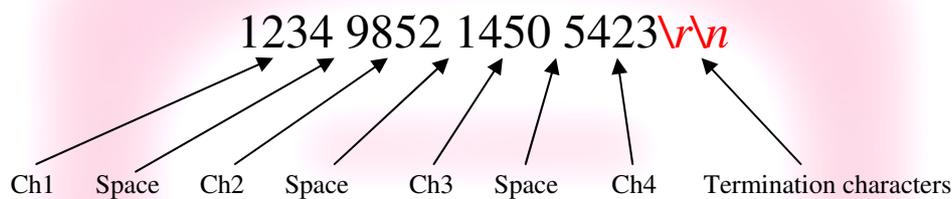

With “BIN ON” the values are transmitted as a continuous binary data stream without any termination or control character. For this reason it is important to keep synchronization with the data flux otherwise it is possible, if some bytes are missed, to corrupt the data and to get false readings from the instrument. If this happens, data acquisition must be stopped, with the “ACQ OFF” command, and then restarted again slowing down the data throughput rate or increasing the reliability of the communication channel. The speed of data transmission is directly given by the selected integration time and the “half mode” selection (see “HLF Command” paragraph for more information).

In binary mode the data output format consists of a continuous sequence of 12 bytes, three bytes for each channel with MSB first:

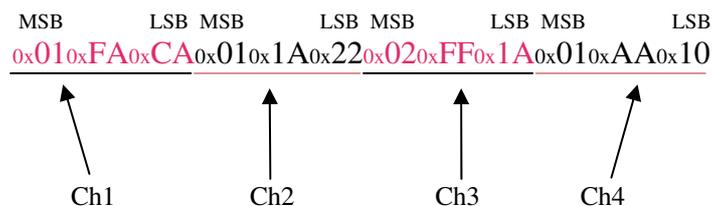

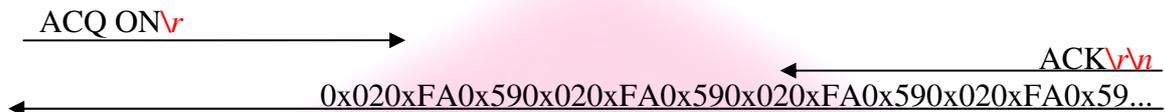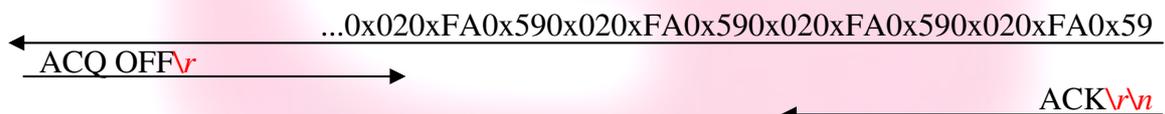Examples:*ACQ ON example with ASCII data:**ACQ ON example with BINARY data:**ACQ OFF example with ASCII data:**ACQ OFF example with BINARY data:**ACQ ? example:*

2.2.2 “BDR” Command

The purpose of the BDR command is to set the baudrate between the AH401B and the communication module. The following table shows the allowed baudrates:

Command	Baudrate
BDR 921600	921600
BDR 460800	460800
BDR 230400	230400
BDR 115200	115200
BDR 57600	57600
BDR 38400	38400
BDR 19200	19200
BDR 9600	9600

By default the instrument boots with a 921600 baudrate setting.

Since the instrument immediately changes the baudrate when the BDR command is accepted, no acknowledgement string (ACK) is sent back to the host. The user should verify successful execution of the command by sending another command after changing the host baudrate (e.g.: “BDR ?”).

The baudrate change procedure strictly depends on the communication interface module used. Please refer to the “Communication Module Interfaces” chapter for more information.

Examples:

BDR example with incorrect parameter:

BDR 960000\r

NAK\r\n

BDR example with correct parameter:

BDR 9600\r

No acknowledgement from the AH401B but the baudrate has been changed. In order to verify that the baudrate was effectively changed the “BDR ?” command should be sent.

BDR ? example:

BDR ?\r

BDR 921600\r\n



2.2.4 “GET” and “?” Command

The purpose of the GET command is to read back a single snapshot of the four channels values. The “?” command is a useful shortcut fully equivalent to the GET command. The format of the returned values depends on the “binary mode” setting (refer to “BIN Command” paragraph).

Examples:

GET ? example with ASCII data:

```
GET ?\r → 8232 43567 9803 7996\r\n ←
```

? example with ASCII data:

```
?\r → 8232 43567 9803 7996\r\n ←
```

GET ? example with binary data:

```
GET ?\r → 0x020xFA0x590x020xFA0x590x020xFA0x590x020xFA0x59 ←
```

? example with binary data:

```
?\r → 0x020xFA0x590x020xFA0x590x020xFA0x590x020xFA0x59 ←
```

2.2.5 “HLF” Command

The purpose of the HLF command is to select whether to process data from both integrator circuits (maximum speed) or only from one integrator circuit (best noise performance) of the AH401B.

In order to explain the functionality of the HLF command it is useful to describe how the integrator input stage operates. A basic feature of the AH401B is that each input channel consists of two parallel integrators (A and B) so that the current-to-voltage integration can be performed continuously in time: when one of the integrators (i.e. A) is digitized by the ADC, the input channel is switched to the other integrator circuit (i.e. B) and the input current can thus be continuously integrated. At the end of the integration time, also the second integrator (i.e. B) is disconnected from the input and the ADC can start the digitization of the integrated current from B. At the same time, the input is connected to the previous integrator (i.e. A) and therefore the current can continue to be integrated by A.

This sequence is continuously repeated as long as the AH401B is in acquisition mode (“ACQ ON”). Dead time in the data throughput is thus avoided and the sampled values are continuously sent to the host. This mode of operation (“HLF OFF” - *default*) is useful when the maximum sampling rate is required since at the end of each integration cycle a digitized data set is sent to the host PC. The drawback is a slightly higher noise level on the sampled values due to the integrator capacitors mismatch between A and B and to the charge injection introduced by the internal switches.

If lowest noise performance is of uttermost importance, the “HLF ON” command must always be used. In this operation mode only the integrated current of one integrator (i.e. A) is sampled, digitized and sent to the host. Using only and always the same integrator any capacitors mismatch and charge injection is avoided and the best signal to noise ratio is achieved. The operating sequence of this mode is as follow: the current is integrated on integrator A for the integration time (i.e. 100ms), then the value is digitized and sent to the host. During the following integration time (i.e. 100ms) no data is digitized (the value on the integrator B is discarded) and then the sequence repeats itself. Therefore a set of data is sent to the host every two integration times (i.e. 200ms). The drawback of this mode is that only “half” sampled values are sent to the host and hence the sampling rate is halved.


Please note that the data rate throughput is directly related to the integration time and the “half mode” selection. For example, setting the integration time to 10ms and the “half mode” to disabled (HLF OFF) generates a data stream at 100Hz. Whereas, setting the integration time to 10ms and the “half mode” to enabled (HLF ON), generates a data stream at 50Hz.




Examples:

HLF ON example:

HLF ON\r

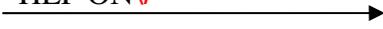

ACK\r\n

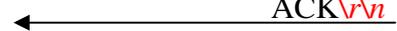

HLF ? example:

HLF ?\r



HLF ON\r\n



2.2.6 "ITM" Command

The purpose of the ITM command is to set the integration time of the AH401B. The parameter field is time expressed in hundreds of microseconds, i.e. if a 100ms integration time is needed the parameter must be set to 1000. The range of the integration time spans from 1ms to 1s, so the parameter range goes from 10 to 10000. At power-up the AH401B integration time is set to 100ms (ITM 1000 - *default*) by default.

Examples:

ITM example for 1ms integration time:

ITM 10\r → ← ACK\r\n

ITM example for 200ms integration time:

ITM 2000\r → ← ACK\r\n

ITM ? example:

ITM ?\r → ← ITM 1000\r\n

2.2.7 “NAQ” Command

The purpose of the NAQ command is to read a fixed number of samples, ranging from 1 to 20000000, without having to manually stop the acquisition (i.e. “ACQ OFF” command) when the desired number of samples has been read. This function has to be used in combination with the ACQ command (refer to “ACQ Command” paragraph) or the TRG command (refer to “TRG Command” paragraph) in order to start the acquisition. The format of the returned values depends on the “binary mode” setting (refer to “BIN Command” paragraph). The NAQ function is disabled when the number of samples is set to 0.

Examples:

NAQ example for 4 samples:

NAQ 4\r → ← ACK\r\n

ACQ ON example with ASCII data:

ACQ ON\r → ← ACK\r\n
 ← 3124 12001 1222 10\r\n
 ← 3143 11991 1342 20\r\n
 ← 3185 12891 1349 25\r\n
 ← 3213 10091 1448 35\r\n

NAQ ? example:

NAQ ?\r → ← NAQ 4\r\n

2.2.8 “RNG” Command

The purpose of the RNG command is to set the full scale range of the AH401B. The parameter field is a number representing the selected charge integration range as reported in the following table:

	Full Scale	Equivalent Current FS @ 1ms	Equivalent Current FS @ 1s
RNG 0	1.8nC	1.8μA	1.8nA
RNG 1	50pC	50nA	50pA
RNG 2	100pC	100nA	100pA
RNG 3	150pC	150nA	150pA
RNG 4	200pC	200nA	200pA
RNG 5	250pC	250nA	250pA
RNG 6	300pC	300nA	300pA
RNG 7	350pC	350nA	350pA

The parameter value can be set from 0 to 7. At power up the AH401B range is set to 1 (*default*), corresponding to a full scale range of 50pC.

Examples:

RNG example for 100pC Full Scale Range:

RNG 2\r → ← ACK\r\n

RNG example for 1.8nC Full Scale Range:

RNG 0\r → ← ACK\r\n

RNG ? example:

RNG ?\r → ← RNG 3\r\n

2.2.9 “SUM” Command

The purpose of the SUM command is to add the values of “N” data samples configured with the NAQ command (refer to “NAQ Command” paragraph) and hence get a single channel value representing the summed samples. This function has to be used in combination with the NAQ command (refer to “NAQ Command” paragraph). The format of the returned values depends on the “binary mode” setting (refer to “BIN Command” paragraph).

Examples:

SUM example with 4 samples configured:

NAQ 4\r


SUM ON\r


ACQ ON example with ASCII data:

ACQ ON\r


SUM ? example:

SUM ?\r


2.2.10 “TRG” Command


The purpose of the TRG command is to enable/disable the “trigger mode” of the AH401B. The “trigger mode”, when enabled (TRG ON), allows synchronizing the picoammeter data acquisition with an external event (for instance laser pulses). The “trigger mode” operation is described in the next paragraph.

As soon as the TRG ON command is received the AH401B stops data acquisition and waits to receive, on the corresponding TRIGGER pin (see “I/O Connectors” chapter), a falling edge signal. As soon as this signal is detected, the AH401B starts to acquire data and immediately and continuously sends it to the host until a new TRIGGER signal is received on the corresponding input. If the NAQ function is enabled (refer to “NAQ Command” paragraph), the acquisition automatically stops after the configured number of samples are transmitted and the instrument waits for a new TRIGGER signal. Moreover, if the NAQ function is used in combination with the SUM mode (refer to “SUM Command” paragraph) only the summed values of the samples are returned. This behavior continues until the “trigger mode” is disabled with the “TRG OFF” (*default*) command (see “TRIGGER Signal” paragraph in the “I/O Connectors” chapter).


The format of the data sent depends on the “binary mode” setting that can be changed using the “BIN” command (see “BIN Command” paragraph).

Examples:


TRG example with ASCII data:

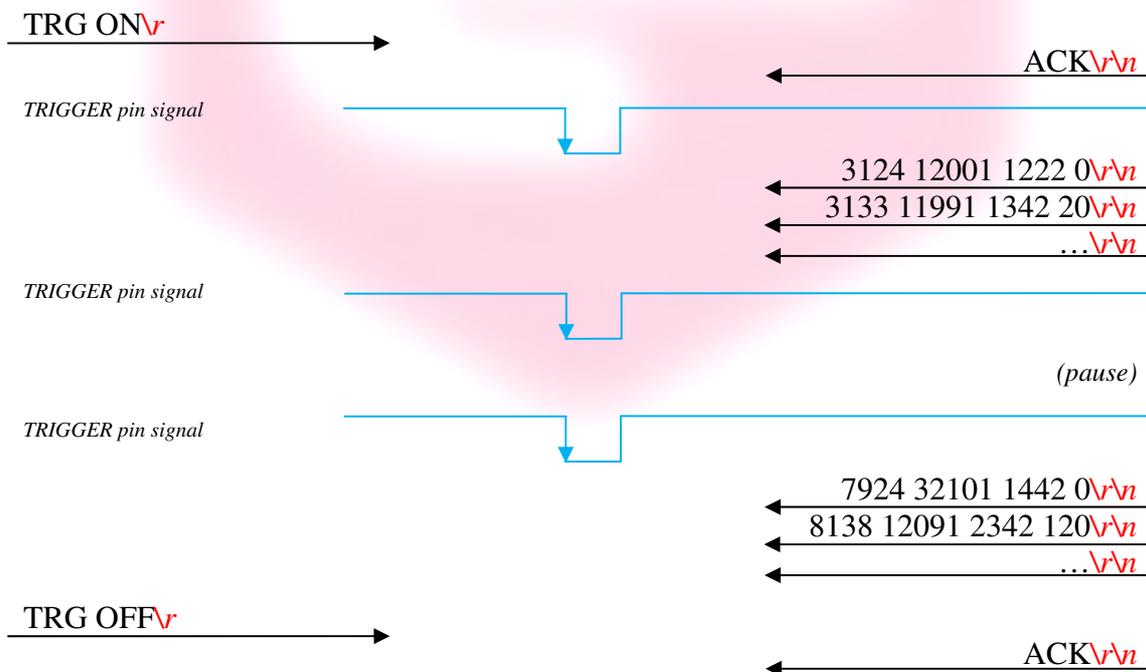


TRG example with NAQ function enabled (ASCII data):

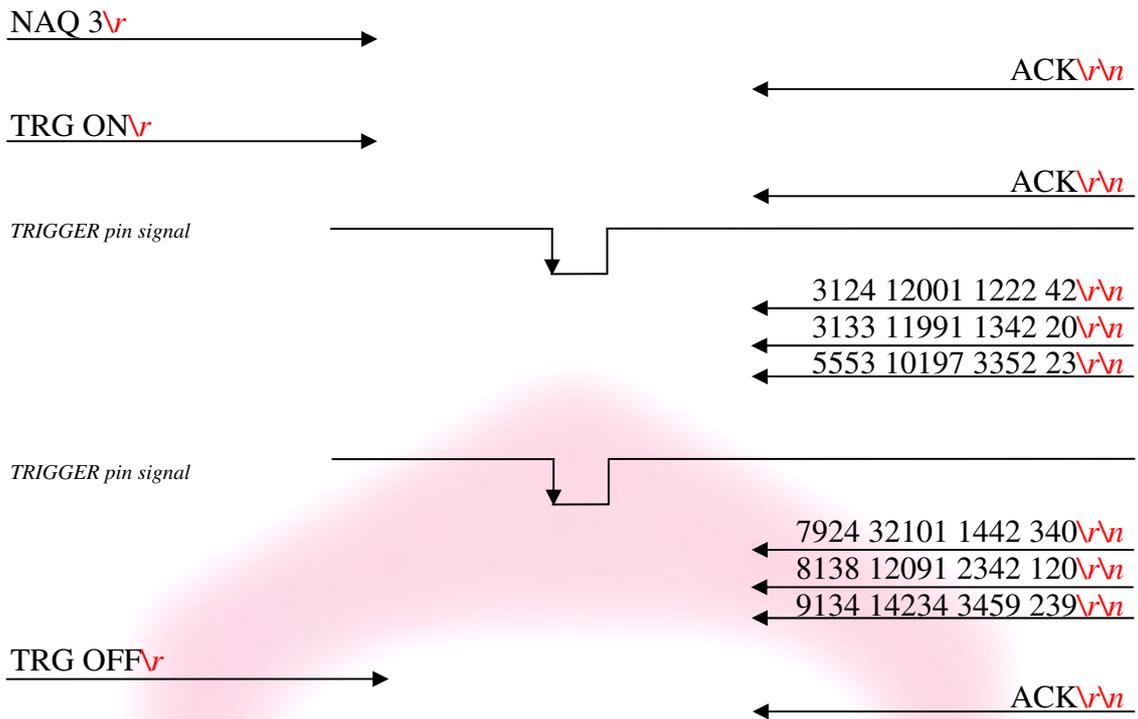

TRG example with NAQ and SUM functions enabled (ASCII data):

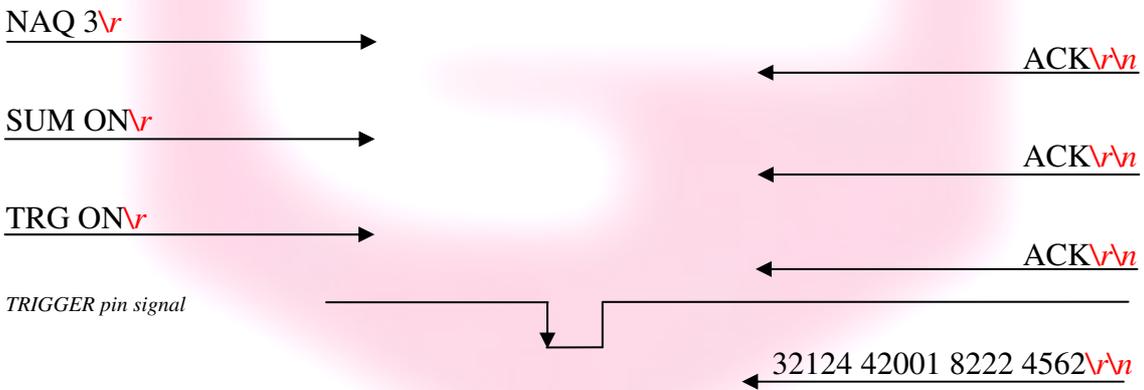

TRG ? example:

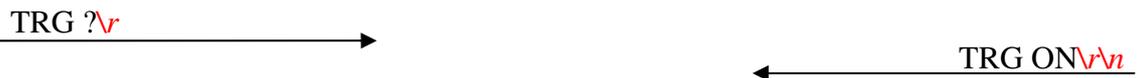

2.2.11 “VER” Command

The purpose of the VER command is to report the firmware version currently installed in the AH401B on-board processor.

Examples:

VER ? example:

VER ?*r*

VER 2.1.0*\r\n*

2.3 Command Table Summary

Command	Purposes	Parameters	Power-Up default value
ACQ	Start continuous acquisition	ON	
	Stop continuous acquisition	OFF	<i>default</i>
	Query acquisition status	?	
BDR	Set baudrate	921600	<i>default</i>
	Set baudrate	460800	
	Set baudrate	230400	
	Set baudrate	115200	
	Set baudrate	57600	
	Set baudrate	38400	
	Set baudrate	19200	
	Set baudrate	9600	
	Query baudrate setting	?	
BIN	Select binary mode	ON	
	Select ASCII mode	OFF	<i>default</i>
	Query "binary mode" status	?	
GET	Read a single snapshot	?	
?			
HLF	Select half mode	ON	
	Select full mode	OFF	<i>default</i>
	Query "acquisition mode" status	?	
ITM	Set integration time	1000	<i>default</i>
	Set integration time	10 to 10000	
	Query integration time	?	
NAQ	Read a fixed number of samples	1 to 20000000	
	Disable NAQ function	0	<i>default</i>
RNG	Set Full Scale Range to 1.8nC	0	
	Set Full Scale Range to 50pC	1	<i>default</i>
	Set Full Scale Range to 100pC	2	
	Set Full Scale Range to 150pC	3	
	Set Full Scale Range to 200pC	4	
	Set Full Scale Range to 250pC	5	
	Set Full Scale Range to 300pC	6	
	Set Full Scale Range to 350pC	7	
	Query Full Scale Range	?	
SUM	Sum the NAQ configured samples	ON	
	Disable Sum mode	OFF	<i>default</i>
TRG	Set trigger mode	ON	
	Disable trigger mode	OFF	<i>default</i>
	Query "trigger mode" status	?	
VER	Query firmware version	?	

2



3. Communication Modules

This chapter describes the currently available communication modules (RS232Piggy, RS422/485Piggy, USBPiggy and ETHERPiggy) and how to install and configure them.

3.1 Module Hardware Installation

The AH401B is shipped with a communication module already installed as shown in Fig. 2.



Fig. 2 - The AH401B with an Ethernet communication module

In order to change the communication module type follow these steps:

- Switch off the instrument;
- Unscrew the module retaining screw;



- Gently pull out the communication module using the retaining screw;



- Insert the new module. Be careful to verify that the module connector and the AH401B receptacle connector are properly aligned;



- Fix the retaining screw;
- Switch on the instrument.

Now the new communication module should be installed and running. To properly set-up each module, please follow the software configuration instructions in the corresponding paragraph.

3.2 “ETHERPiggy” Module

The “ETHERPiggy” module is based on the Lantronix Xport® embedded Ethernet device server. To properly configure the Xport® device the user must configure the IP address, the communication protocol and all other network parameters needed by the AH401B. Please, refer to the Xport® documentation and configuration software available at the Lantronix site www.lantronix.com for a complete description.

3.2.1 “ETHERPiggy” IP Address Assign

Every “ETHERPiggy” module is shipped with a default IP address as indicated on a label on the module itself (see Fig. 3).



Fig. 3 - "ETHERPiggy" Module with label

Before assigning a new IP address it is required to install, on the host PC, the “DeviceInstaller” software that can be downloaded from the Lantronix website www.lantronix.com.


Even if the AH401B can be connected on the global LAN, a point-to-point Ethernet connection is strongly recommended in order to insure maximum speed performance and to avoid possible communication problems. This implies that the host PC and the “ETHERPiggy Module” should reside on the same ethernet subnet, i.e.:

Host PC address: 192.168.10.100 → ETHERPiggy address: 192.168.10.200


Please note that in a point-to-point direct connection a twisted Ethernet cable must be used.

In order to assign the IP address to the “ETHERPiggy” module the next steps must be followed:


- Connect the host PC and the AH401B with a twisted Ethernet cable;
- Switch on the AH401B;
- Verify that the “Link LED” on the RJ45 connector is on (amber for a 10Mbps connection, or green for a 100Mbps connection) and the corresponding “Network Connection” is active on the host PC;

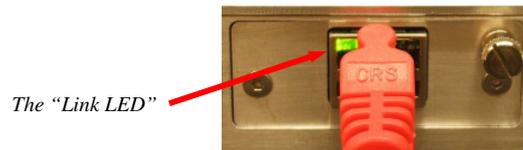


- Launch the “DeviceInstaller” program;
- Select the XPort device where you want to change the IP address;

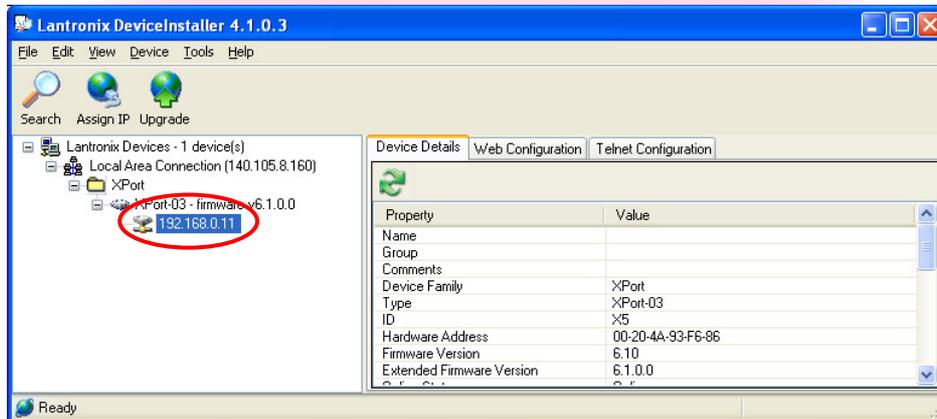


- Click on the “Assign IP” icon;

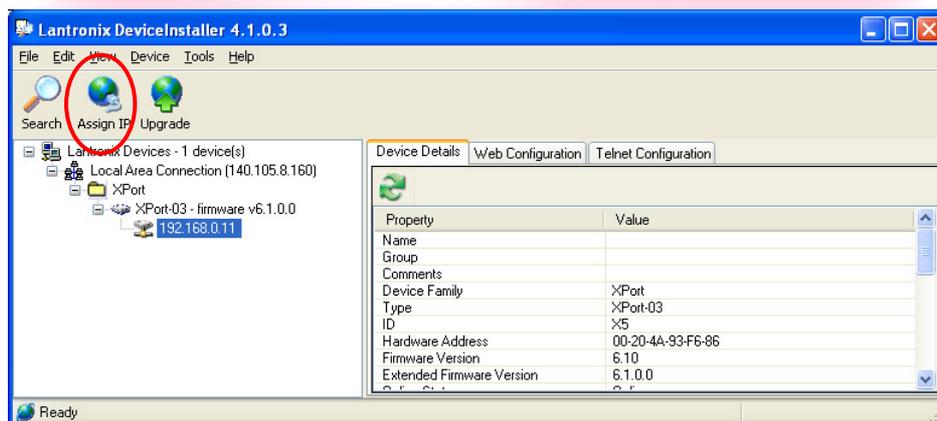


- Select “Assign a specific IP address” and then click “Next”;

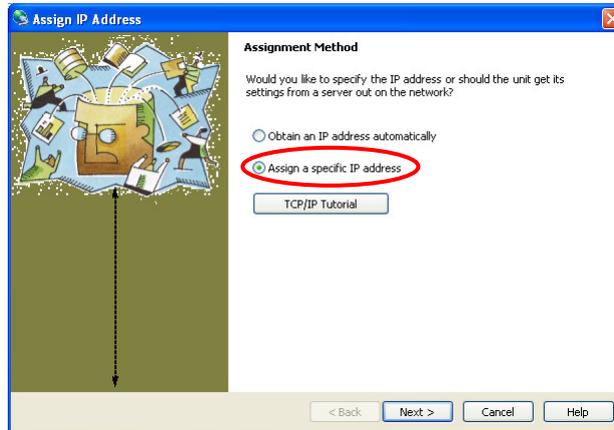

- Set the “IP address”. Click “Next”;

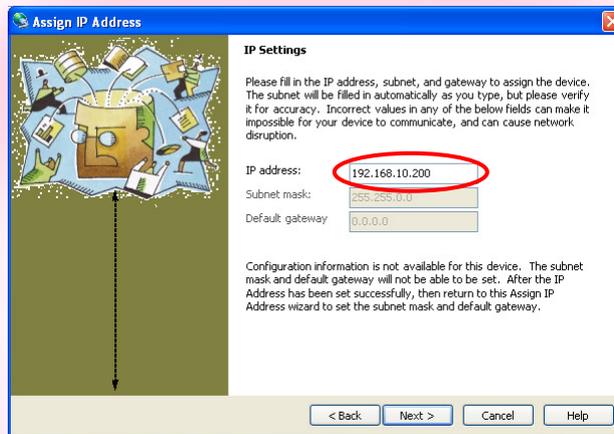

- Click the “Assign” button;

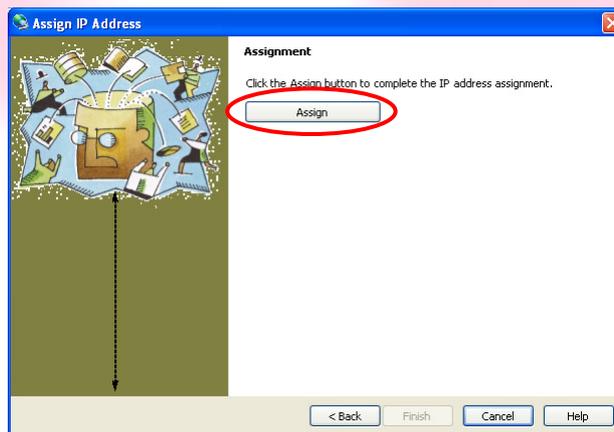

- Wait the “Assignment” then click “Finish”.

Now the new module IP address should be assigned and the success of the operation can be verified on the “DeviceInstaller” window.

3.2.2 “ETHERPiggy” Software Configuration


After assigning the IP address, the user must configure some other parameters on the Xport® device. There are two possibilities to configure such parameters:

- using the web server interface;
- using a telnet connection.

In this manual we describe only the web interface procedure, if you need to configure via the telnet connection please refer to the Xport® user guide.


Follow the next steps to properly setup the ETHERPiggy communication module:

- Select the “Web Configuration” tab, then click the “Go” button. Be sure to disable the “proxy server” option in your “Internet Options” application: go to the “Control Panel” folder, open the “Internet Options”, then click on the “Connections” tab, go to the “LAN settings” and then deselect the “Use a proxy server for your LAN” option;

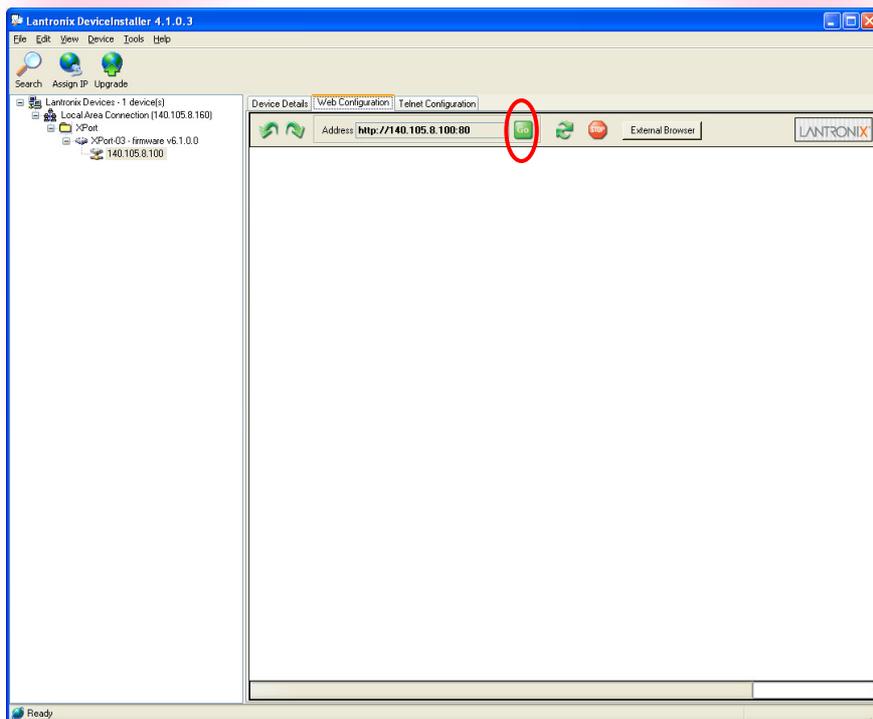


- A dialog window will appear asking for user name and password, simply click the “OK” button;


- Then, the Lantronix web server interface will appear;

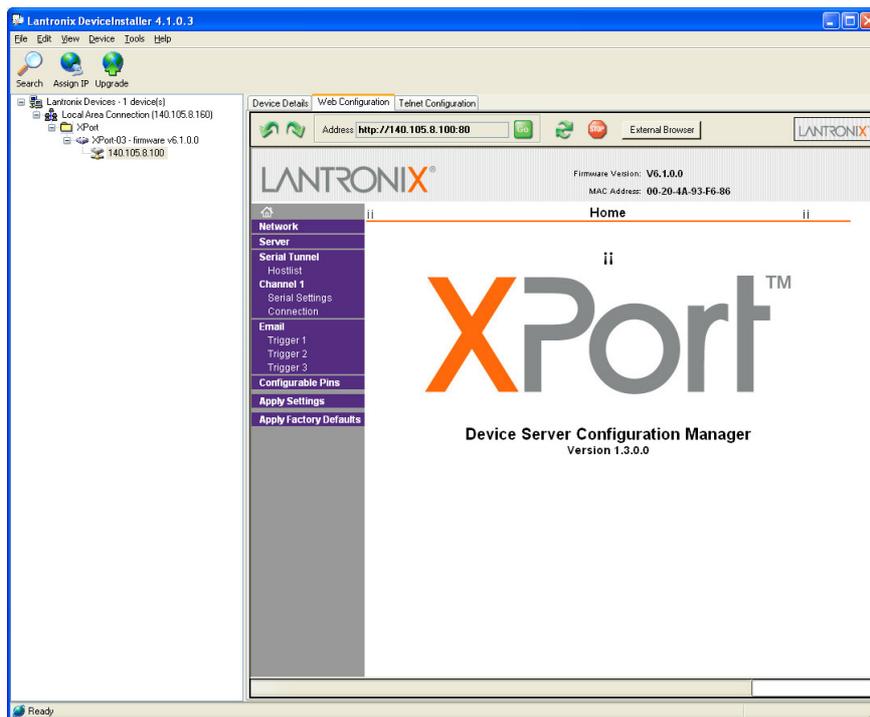


- Click on the “Network” button and set the “Subnet Mask” to “255.255.255.0”; Then click the “OK” button;

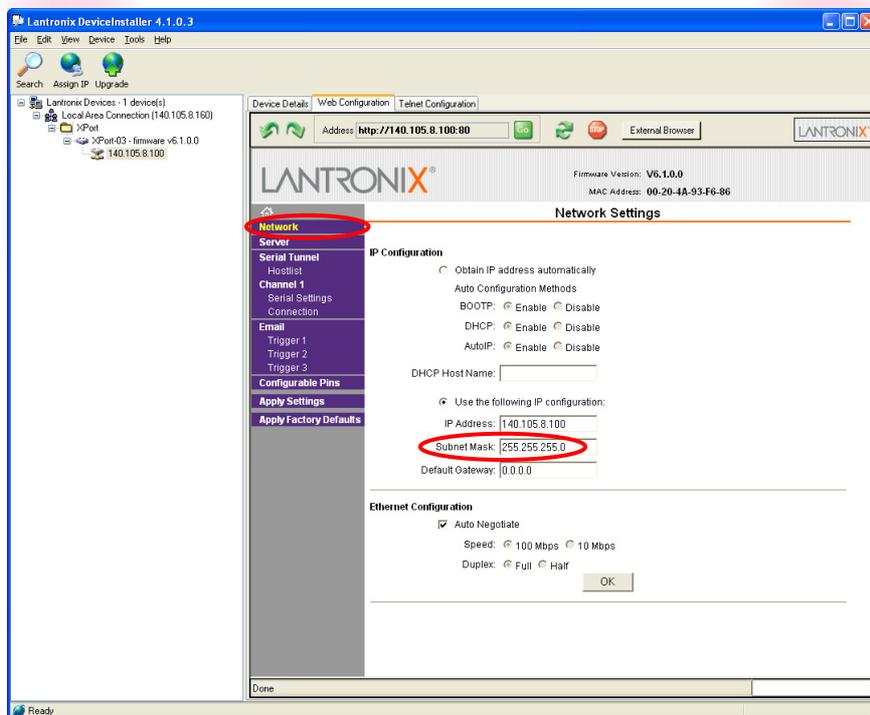


- Click on the “Server” button. Set the “CPU Performance Mode” to “High”. Click the “OK” button;

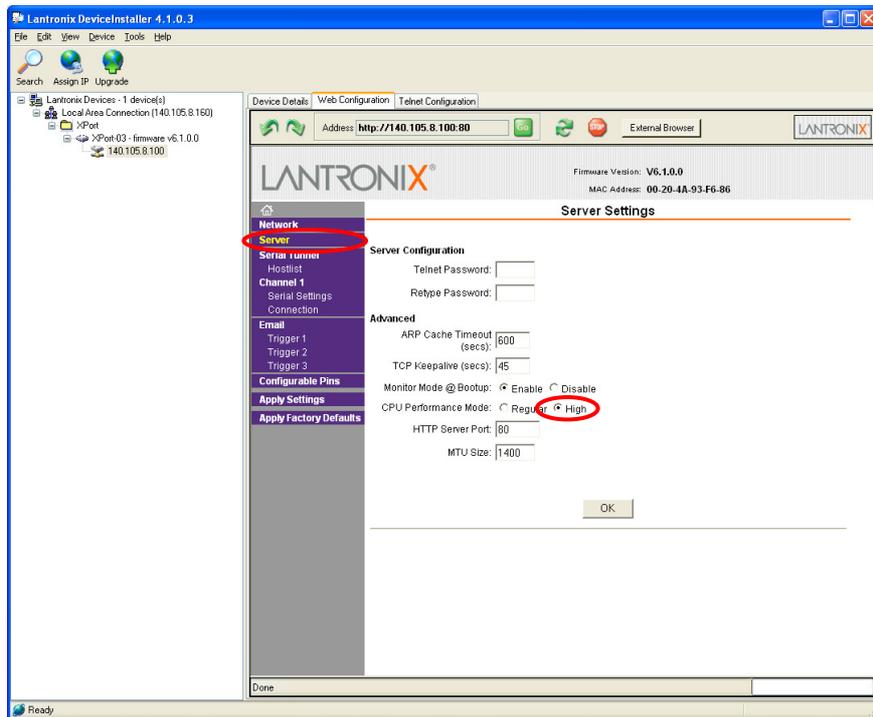

- Click on the “Serial Settings” button. Set the “Baud Rate” to “921600” and the “Flow Control” to “None”. Click the “OK” button;

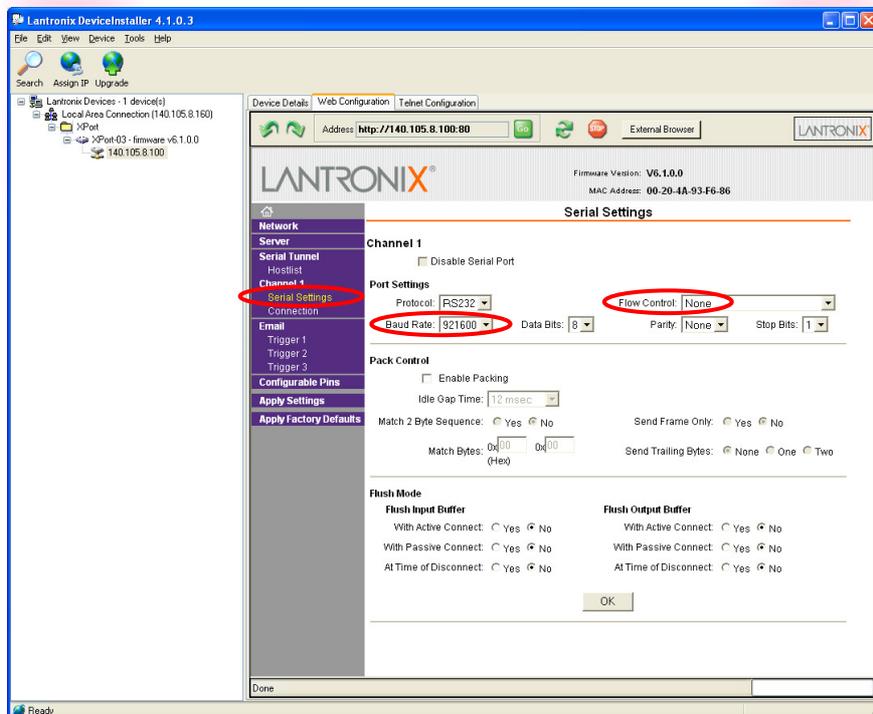

- Click on the “Connection” button. Set the “Protocol” to “TCP” and the “Local Port” to “10001”, this value can be changed as needed. Click the “OK” button;

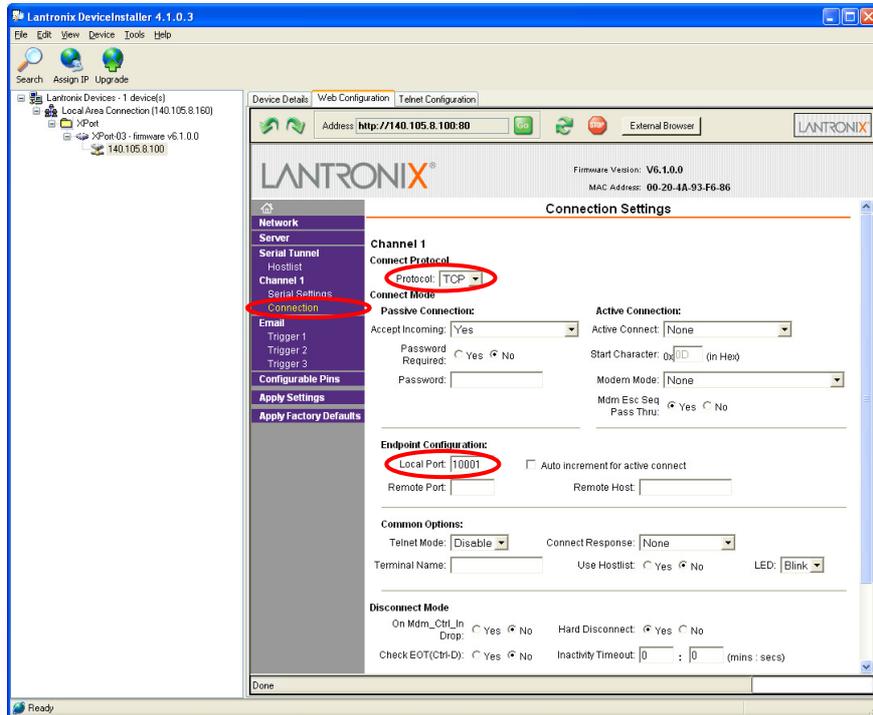

- Click the “Apply Settings” button.

Now the Xport® Ethernet device should be correctly configured to work with the AH401B. If you need to change any other XPort® parameter, please, refer to the Xport® user's guide.

4. I/O Connectors

This chapter describes the I/O connectors, their corresponding pinout and the purposes of each signal.

4.1 Power Connector

In Fig. 4 the input Power Connector is shown with its corresponding pinout. The input voltage range from 9V up to 15V with a max input current of 350mA (depending on the communication module).

The connector type is a LEMO – “EGG.0B.302” series. The corresponding matching connector plug is a LEMO – “FGG.0B.302” series.

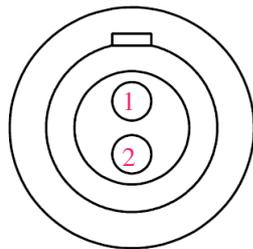


Fig. 4 - Power Connector

Pin	Signal
1	+V _{CC}
2	GND

4.2 I/O Interface Connector

In Fig. 5 the I/O Interface Connector is shown with its corresponding pinout. The signal levels are 5V TTL and CMOS compatible. The minimum output current is 24mA.

The connector type is a DSub 9-pin socket.

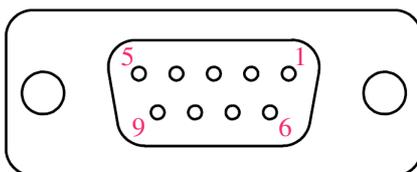


Fig. 5 - I/O Interface Connector

Pin	Signal
1	GND
3	CONV
5	TRIGGER

4.2.1 The CONV Signal

The CONV signal is a copy of the internal Integration Time signal. For example with an Integration Time of 100ms the CONV output will be a square wave with the half-period equal to 100ms. It can be used as a monitor of the internal timing or as synchronization signal for external devices.

The output voltage level is 5V TTL and CMOS compatible, with a minimum output current of 24mA.

4.2.2 The TRIGGER Signal

The purpose of the TRIGGER signal is to synchronize the acquisition of the AH401B with external events (e.g. laser pulses). It works in combination with the TRG command (please refer to “TRG Command” paragraph for more details).

When the “trigger mode” is enabled by software, a falling edge signal on the TRIGGER input starts the acquisition of the picoammeter. As soon as a new TRIGGER signal is received on the corresponding input, the acquisition stops and no data is sent to the host. This behavior continues until the “trigger mode” is disabled by software.

The input voltage level is 5V TTL and CMOS compatible.

5. Technical Specifications

Input Channels	4	
Current Measuring Range	from 50 pA to 1.8 μ A	
Current Polarity	positive only	
Integration Time	from 1ms to 1s	
Data Transfer	up to 1 ksamples/s	
Resolution Bits	20	
Noise (@1ms, 200pC_{FS})	<7ppm	
Communication Modules	RS-232/422/485, USB, Ethernet TCP/IP and UDP	
I/O Signal	CONV output - TRIGGER input	
Supply Voltage	from 9 V to 15 V	
Max Supply Current	RS-232/422/485	100 mA
	USB	150 mA
	Ethernet (TCP/IP and UDP)	350 mA
Dimensions	140 × 110 × 28 mm	
Weight	420 g	
Input Connectors	BNC	